

HEATING MANTLES

Heating Mantle. Simple design and suitable for use in schools universities and laboratories.

Features

- Flameless Heat
- Large contact area to glassware
- Even heat distribution
- Rapid continuous use
- Powder coated for chemical resistance
- Wired in Australia by IEC Pty Ltd



Instructions:

- Power ON and lamp will indicate
- Rotate knob clockwise to increase heating
- An indicator lamp will cycle with the heating element

Temperature: 360 deg c. Surface of heating element.

Note: The first use will cause smoke and create odour for up to an hour until the coating on the element is dissipated. First operation should occur in a well ventilated area.

SIZE	POWER	CODE NO
100ML	220/240 V.AC 50/60Hz 0.5A 100W	GE3277-01
250ML	220/240 V.AC 50/60Hz 0.7A 150W	GE3278-01
500ML	220/240 V.AC 50/60Hz 1.2A 280W	GE3279-01
1000ML	220/240 V.AC 50/60Hz 1.7A 400W	GE3280-01